
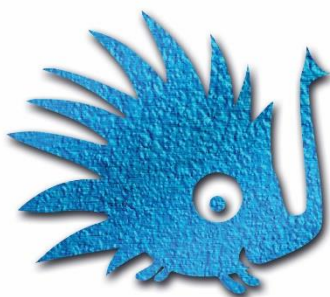


Mouse Wrap

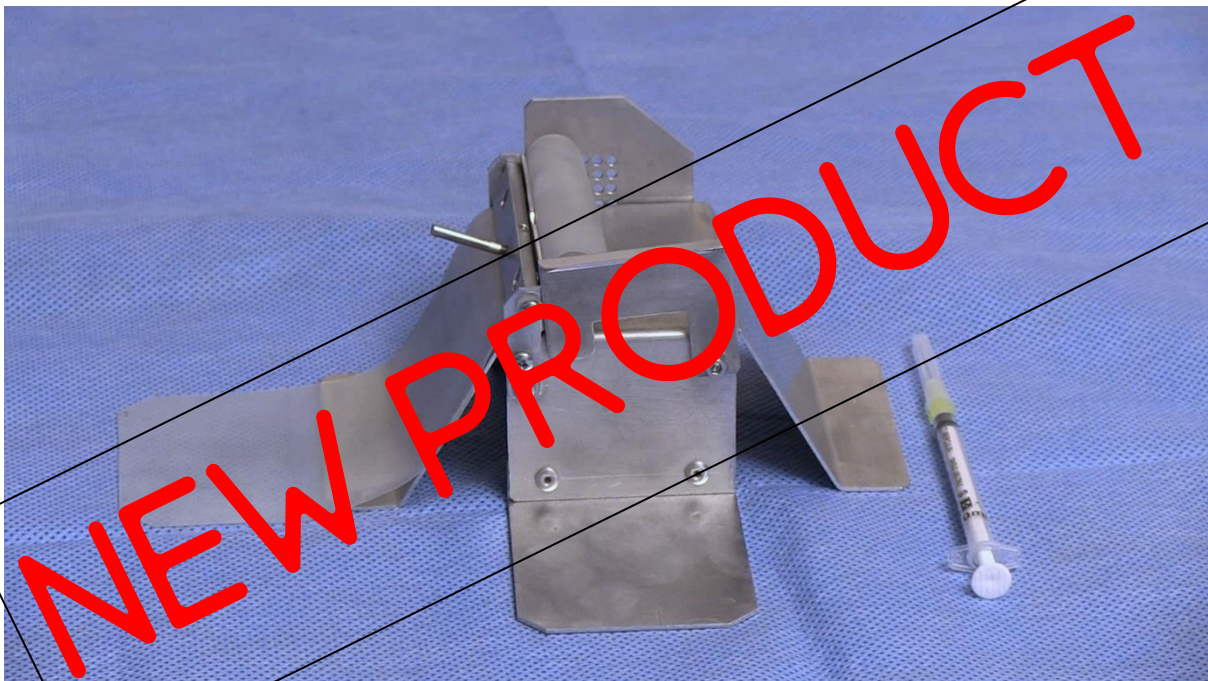

OTAWOG

Rodent Anaesthesia Systems

How to Use the Mouse wrap

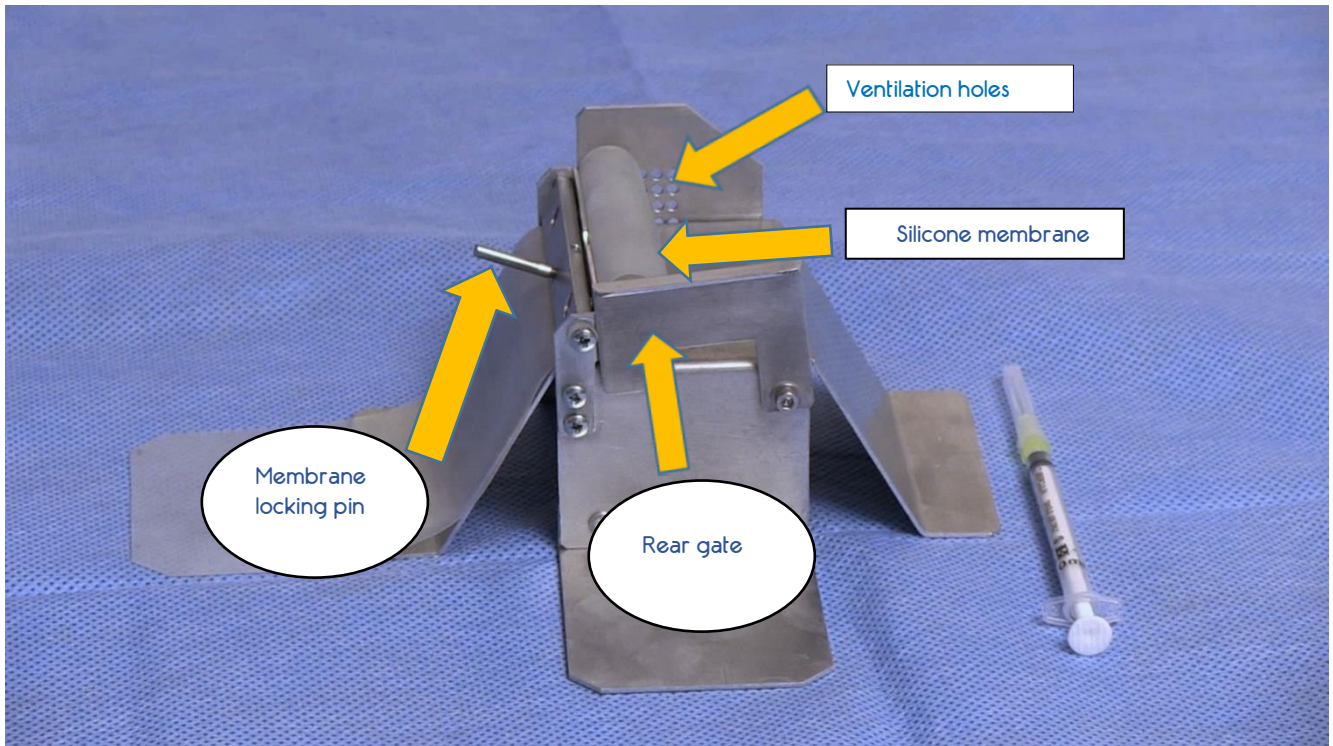


Figure 1: Front view of Mouse Wrap

To wrap a mouse:

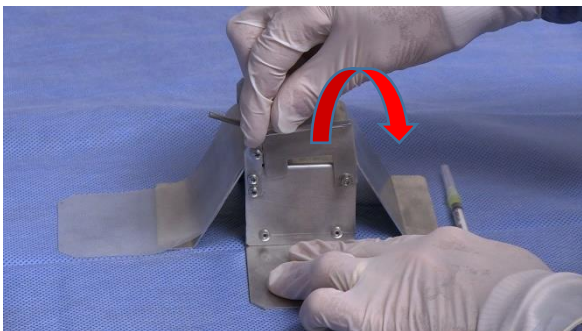


Figure 2: Unlock the rear gate

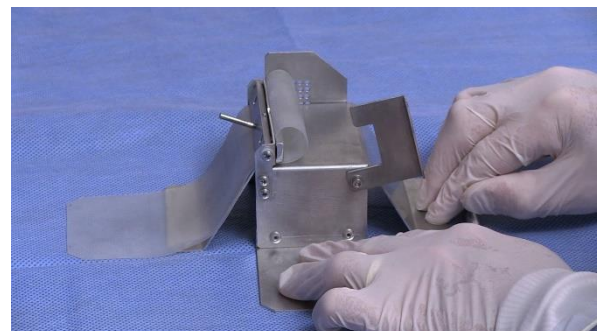


Figure 3: Rear gate unlocked. Shows the silicone tunnel.

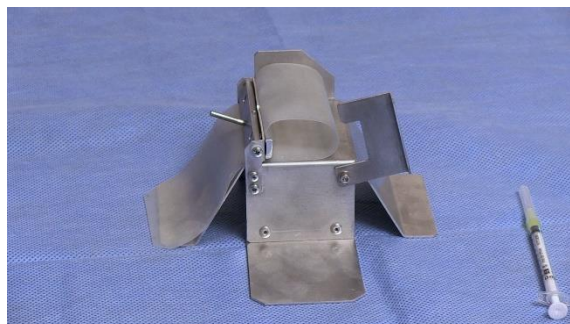


Figure 4: Silicone tunnel has been enlarged so that a mouse can be introduced.

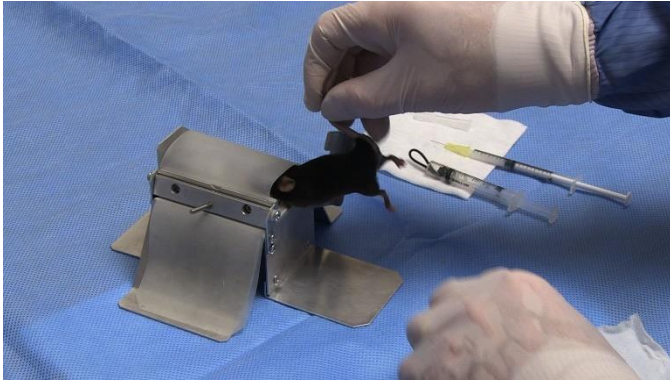


Figure 5: The mouse is shown the tunnel. It will run in

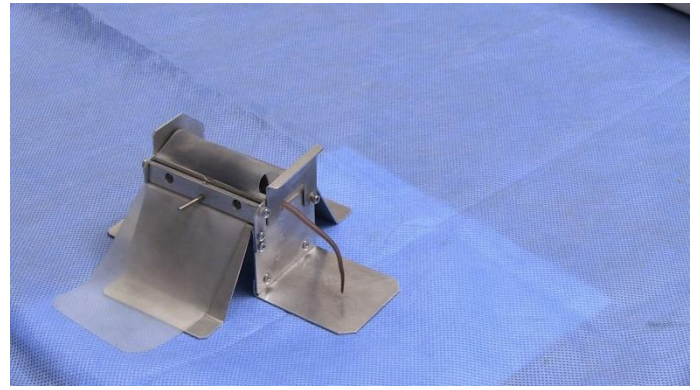


Figure 6: Pull the membrane to snug around the mouse. Lock the tab and replace the rear gate over the tail



Figure 7: The tail is now accessible for IV, tail tip, tattoo or blood collection

To release the mouse



Figure 8: To release the mouse, unlock the tail gate, Unlock the membrane and allow mouse to exit the front or back of the device

Disinfection or sterilisation recommendations:

1. Rinse the Wrap in warm soapy water to remove any gross contamination.
2. The Wrap can then be sprayed with 70% alcohol and left to dry.
3. The Wrap can also be autoclaved (121C for 15 minutes).



www.otawog.com

Dunedin, New Zealand
info@otawog.com